

TENAK LabBird blink patterns

LabBird

Online

Has IP address and TENAK MyLog cloud connection.

- Constant green LED.

Offline

Has either no IP address or no TENAK MyLog cloud connection.

- Constant red LED.

Parring mode

Entered by a short button press.

- Blinking green for 60 s. or until pairing has completed.

Radio region missing

- Alternating blinking red and green.

Firmware update pending

- Alternating blue and green blinks. Scheduled LabLogger firmware upgrade. LabLogger will be updated on next contact with LabBird.

Firmware update in progress

- Constant blue. With normal radio coverage this will take approximately 90 s.

Slave parring

5 s. button push until LED turns blue.

- Blue LED blinks for 60 s. or until paring was completed.

You can always contact us at info@tenak.com with questions, and we will do our best to help you out as fast as possible.

