

TENAK LabLogger blink patterns

LabLogger

Pairing mode

Entered by a short button press, when not already paired.

- Green blinking LED for 60 s.
- On succes, 2 short green blinks and LED turns off.

Sending measurement

- 2 or 3 green blinks (last blink is optional and is ACK from the LabBird).

Manual measurement

Initiated by short button press, on a paired LabLogger.

- 4 green blinks with probe attached (last blink is optional and is ACK from the LabBird).
- 3 green blinks without probe (last blink is optional and is ACK from the LabBird).

Firmware update

- Constant blue light on the LabLogger with serial greater than 20-02-xx-xx.
- No LED indication on other LabLoggers.

Reset

Entered by pressing button for 10 s.

- First no light, then green LED for 5 s. and then it turns red, to indicate button should be released.
- If button is kept pushed LED turns white and reset is canceled.

You can always contact us at info@tenak.com with questions, and we will do our best to help you out as fast as possible.

