

TENAK LabBird

Master / Slave Set-up Guide

Step 1: **Add first LabBird (master) to your account**
You may need the LabBird serial number from the label on its backside. Usually though the serial number can be automatically inferred and filled in for you.

Step 2: **Pair master and slave LabBirds**
Press and hold the button on the LabBird that will be operating as a slave. The button should be pressed for approx. 5 sec. until the LED turns blue and then release button. The LED will now be flashing blue.

Now short press the button on the master LabBird. After a couple of seconds the flashing LED on both LabBirds should stop and the LabBirds are now in master/slave mode.

Confirm master/slave mode on the device-list at mylog.tenak.com where the slave LabBird should be grouped with the master and have a blue slave-label next to the name

If more slave LabBirds are to be added to the setup repeat this step for the next slave LabBird.

Step 3: **Add LabLoggers to master LabBird**
LabLoggers has to be paired to the master with the normal procedure.

You can also contact us at info@tenak.com, and we will do our best to help you out as fast as possible.

